

# 梅特勒-托利多

## PL602-L 电子天平简易操作指南

METTLER TOLEDO

PL602-L 称量参数：

最大量程：610g

可读性：0.01g

秤盘尺寸：直径 120mm

校正砝码：500g F1







操作天平之前的准备工作：






- 1、 选择稳定无振动水平的工作台位置摆放天平。
- 2、 通过天平的调节水平脚，使电子天平水平泡在中央位置。
- 3、 通电预热，精密天平一般 30 分钟。
- 4、 选择 500g 砝码对天平校正。
- 5、 校正完毕，天平就可以称量了。

操作键功能一览

该系列天平具有两种操作方式:称量模式和菜单模式。根据所选择的操作方式和按键时间的长短，各键有不同的含义。



称量模式下的操作键功能	
短时间按键 	长时间按键 
<b>1/10d</b> <ul style="list-style-type: none"><li>实际分度值减小</li></ul>	<b>Cal</b> <ul style="list-style-type: none"><li>调整 (校准)</li></ul>
<b>On</b> <ul style="list-style-type: none"><li>开机</li><li>清零 / 去皮</li><li>取消功能</li></ul>	<b>Off</b> <ul style="list-style-type: none"><li>关机</li></ul>
 <ul style="list-style-type: none"><li>转换</li><li>改变设置</li></ul>	<b>F</b> <ul style="list-style-type: none"><li>功能调用 所需功能须在菜单中激活，否则在显示屏上将出现“F nonE”</li></ul>
 <ul style="list-style-type: none"><li>通过接口传输称量数据到激活的打印机</li><li>数据设置确认</li></ul>	<b>Menu</b> <ul style="list-style-type: none"><li>菜单调用 (按住键不放，直到出现“MENU”)</li></ul>

菜单模式下的操作键功能	
短时间按键 	长时间按键 
<b>1/10d</b> <ul style="list-style-type: none"><li>改变设置</li><li>显示数值减小1位</li></ul>	<b>1/10d</b> <ul style="list-style-type: none"><li>数值快速减小</li></ul>
<b>C</b> <ul style="list-style-type: none"><li>退出菜单 (不保存退出)</li></ul>	-
 <ul style="list-style-type: none"><li>改变设置</li><li>显示数值增加1位</li></ul>	 <ul style="list-style-type: none"><li>数值快速增加</li></ul>
 <ul style="list-style-type: none"><li>选择下一个菜单项</li></ul>	<b>Menu</b> <ul style="list-style-type: none"><li>保存并且退出菜单</li></ul>

# 快速使用指南




点击键



按住不放，直到出现所需提示

▼ 自动步骤

## 开机

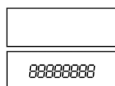

0.00 g

## 关机



OFF



## 基础称量



0.00 g

182.03 g

250.00 g

## 外部调整（校准）



CAL



Cal 2000.00



Cal ---



Cal 0.00



Cal ---



CAL done



0.00 g

## 称量值回忆\*



0.00 g

26.72 g



0.00 g

\* 26.72 g



0.00 g

## 计件\*



0.00 g



12.07 g



PCS 10



PCS 20



20 PCS



244 PCS



147.25 g

## 百分比称量\*



0.00 g

12.07 g



SEE 100%



100.00 %



10.160 %

## 去皮



0.00 g

12.101 g



0.00 g



95.97 g



- 12.101 g

## 称量单位转换\*




64.71 g

647.10 mg

菜单

总述

在菜单中，您可以改变称量单位，选择其他的功能以及进行各种不同的设置。您可以在第 4.3 节中找到各菜单项的说明。

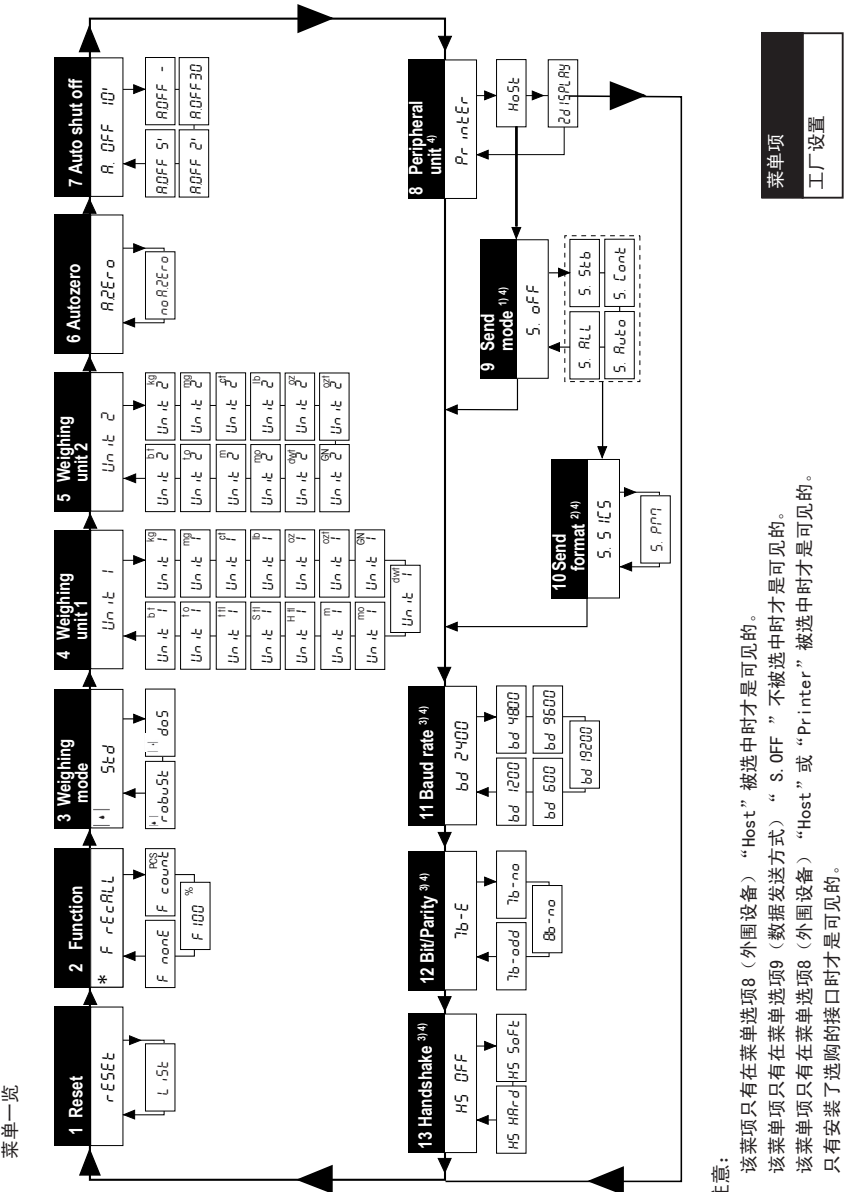

- 注意:
- 1) 该菜单只有在菜单选项8（外围设备）“Host”被选中时才是可见的。
  - 2) 该菜单项只有在菜单选项9（数据发送方式）“S.OFF”不被选中时才是可见的。
  - 3) 该菜单项只有在菜单选项8（外围设备）“Host”或“Printer”被选中时才是可见的。
  - 4) 只有安装了选购的接口时才是可见的。

梅特勒-托利多国际贸易(上海)有限公司  
上海市桂平路 589 号 邮编: 200233  
电话: 021-64850435 传真: 021-64853351  
服务热线: 4008 878 989  
<http://www.mtchina.com> E-mail: [ad@mt.com](mailto:ad@mt.com)

**METTLER TOLEDO**

The logo graphic consists of a series of parallel, slightly curved lines in a light green color, forming a triangular shape that points downwards and to the right. The lines are evenly spaced and have a slight gradient, giving it a sense of depth and movement.